10th World Meeting on Pharmaceutics Biopharmaceutics and Pharmaceutical Technology | Glasgow 4th – 7th April 2016 www.worldmeeting.org



STUDYING THE IMPACT OF LUBRICATION ON TABLET PROPERTIES USING STYL'ONE SUSANA NIETO-BOBADILLA

BRUNO LECLERCQ

THIERRY MENARD



INTRODUCTION

Lubrication plays an important role during tablet manufacturing, especially during ejection and take-off. It reduces the friction between the tablet and the metal surface of the die and punches avoiding common problems as capping, lamination and adhesion to the tooling [1].

However, it is known that lubricants can have negative effects on tablet physical properties; they can interfere with the bonding forces of the particles to be compressed, and might also hinder water penetration, resulting in defects on tablet mechanical strength and dissolution properties [2,3].

The selection of the lubricant and its concentration, as well as the type, speed and time of blending need to be studied during tablet formulation. In the same way, the impact of the press feeding system on the powder lubrication must be evaluated early in R&D. All these parameters affect the lubrication process and consequently the tablet properties. In this context, Styl'One[™], a uniaxial tablet press, permits to rapidly evaluate the impact of lubrication on tablet mechanical properties using only small quantities of powder.

The purpose of this work was to study the influence of the feeding system type on the tensile strength and the ejection force of different types of formulations lubricated with magnesium stearate (MgSt).

EXPERIMENTAL METHODS





Tablets were made either with microcrystalline cellulose (MCC); lactose or a blend of MCC: lactose (30:70), all lubricated with 0.5% MgSt. Tablets were produced using a Styl'One[™] (MedelPharm) (Fig. 1), equipped with Euro B 11.28 mm flat punches.

Two types of feeding systems were used: a gravity and forced feeding shoe. The speed of the last one was set up at 10, 50 and 100% (24.5; 122.5 and 245 RPM). A study of the compression force was performed on Styl'One[™] using the Analis[™] software (Fig. 2).



Fig. 2: Analis[™] acquisition software





Fig. 1: Styl'One™ tablet press and feeding shoes



RESULTS AND DISCUSSION

The impact of the feeding type and forced feeding speed is less critical for lactose (brittle material) (Fig. 3) than for MCC (plastic material) tablets. Over-lubrication at high speed forced feeding causes a remarkable decrease on tensile strength of MCC tablets and the appearance of capping (Fig. 4). The use of external lubrication (punches & die) can avoid the tensile strength drop (Fig. 5). The ejection force can be used as an indicator of the tablet lubrication state (Fig. 6). If lubrication is not early optimized, severe production problems as powder sticking on punches & die, can occur.



Fig. 6: Ejection forces of tablets made of MCC; lactose or a blend of MCC:lactose (30:70) at different feeding shoe conditions

CONCLUSION

ent feeding shoe conditions

Inadequate lubrication affect several tablet parameters. Over-lubrication has a negative impact on tablet tensile strength that could consequently lead to defects on tablet disintegration and dissolution.

feeding shoe conditions

Hence, it is important to choose the right type of excipient (brittle or plastic) and to consider the potential impact of the feeding type (gravitary or forced feeding) on tablet lubrication before to start a production process.

These parameters can be easily and quickly evaluated using a Styl'One[™] compaction simulator during early R&D formulation to ensure appropriate tablet properties.



tion and internal lubrication (0.5% MgSt) at feeding shoe conditions

REFERENCES

1. Wang, J.; Wen, H. and Desai, D. Lubrication in tablet formulations, Eur. J. Pharm. Biopharm. 75, 1-15 (2010).

2. Kikuta, J. and Kitamori, N. Effect of mixing time on the lubricating properties of magnesium stearate and the final characteristics of the compressed tablets, Drug Dev. Ind. Pharm. 20, 343-355 (1994).

3. Otsuka, M.; Yamane, I. and Matsuda, Y. Effects of lubricant mixing on compression properties of various kinds of direct compression excipients and physical properties of the tablets, Adv. Powder Technol. 15, 477-493 (2004).

MEDELPHARM S.A.S

12, rue des petites Combes, Z.I. Nord • F-01700 Beynost • France